



# The Performance Characteristics of a Newly Available ICP Time-of-Flight Mass Spectrometer

Robert Peile, Dmytro R. Bandura, Lindsay Moore

GBC Scientific Equipment Pty Ltd

12 Monterey Road, Dandenong, Victoria 3175, Australia

Greg Bercovy, GBC Scientific Equipment Inc.

3930 Ventura Drive, Arlington Heights, IL, 60004, USA

**GBC**

OptiMass 8000  
ICP Time-of-Flight Mass Spectrometer

1

## **Objectives**

**To evaluate the performance characteristics of a newly-available commercial ICP Time-of-Flight MS, the GBC OptiMass 8000, in terms of :**

- **Resolution**
- **Dynamic Range**
- **Detection Limits**
- **Effectiveness of Ion Blanking**
- **Mass Response**
- **Oxide and Doubly Charged Species Ratios**
- **Ratio Precision**

2

## GBC OptiMass 8000 ICP-TOFMS



# 3

## Description of the OptiMass 8000

**Solid state 27.12 MHz RF generator.**


**Conventional three cone, water cooled interface with differential pumping.**

**Orthogonal acceleration time-of-flight mass spectrometer.**

- **30,000 full spectra per second ion extraction speed.**
- **SMARTGATE ion blanker.**
- **Discrete dynode multiplier detector.**
- **400 MHz detection system sampling rate.**
- **Dual TDC/TR detection system.**

4

# Schematic of the OptiMass 8000

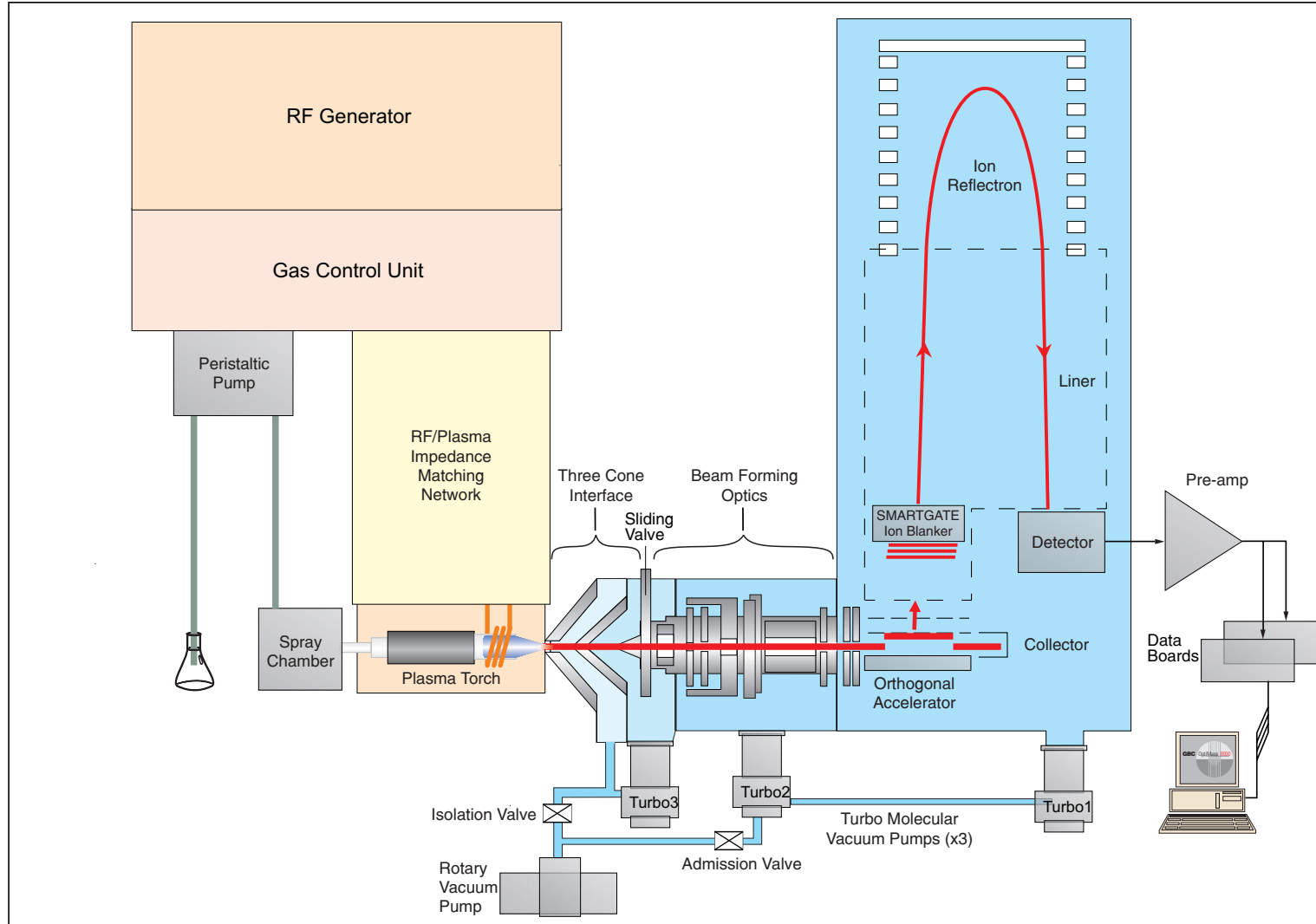

# 5

## Evaluating Resolution

- Resolution in TOF MS is defined as

$$R = M/\Delta M = T/(2\Delta T),$$



where T is time of flight.

As  $T = \text{Const.} \times \sqrt{M}$ , resolution depends on mass

- Resolution of OPTIMASS was evaluated for different masses by 7 x 18 s TDC measurements (integral of 557,056 OA extractions for each measurement) of 10 ppb multi-element RO water solution


# 6

## Typical Resolution

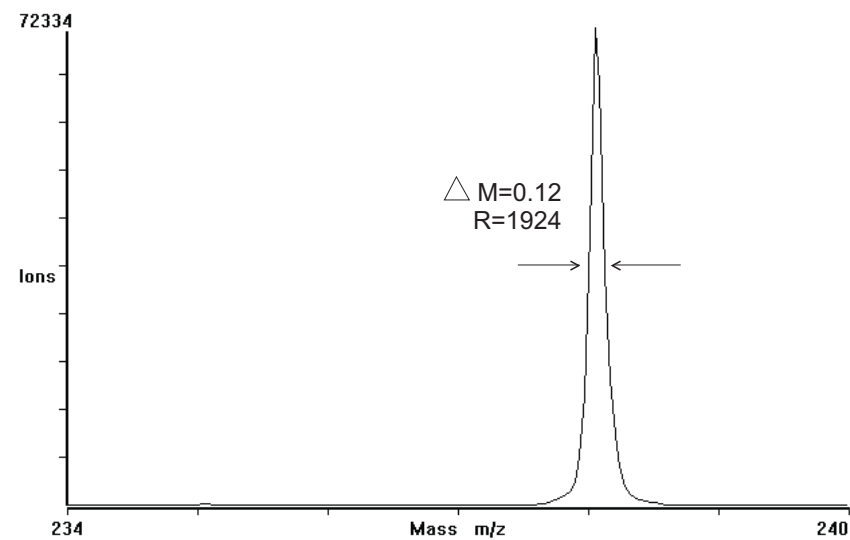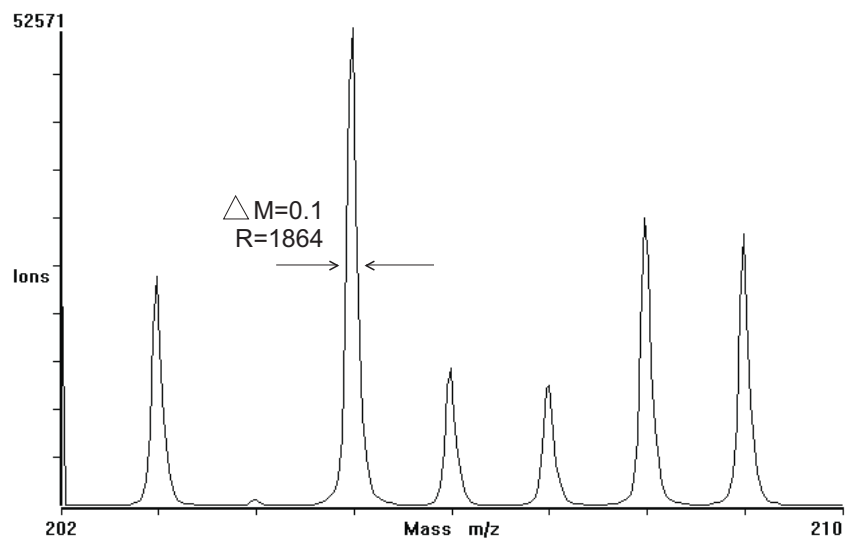

# Resolution vs. mass

Resolution (f.w.h.m.) Versus mass

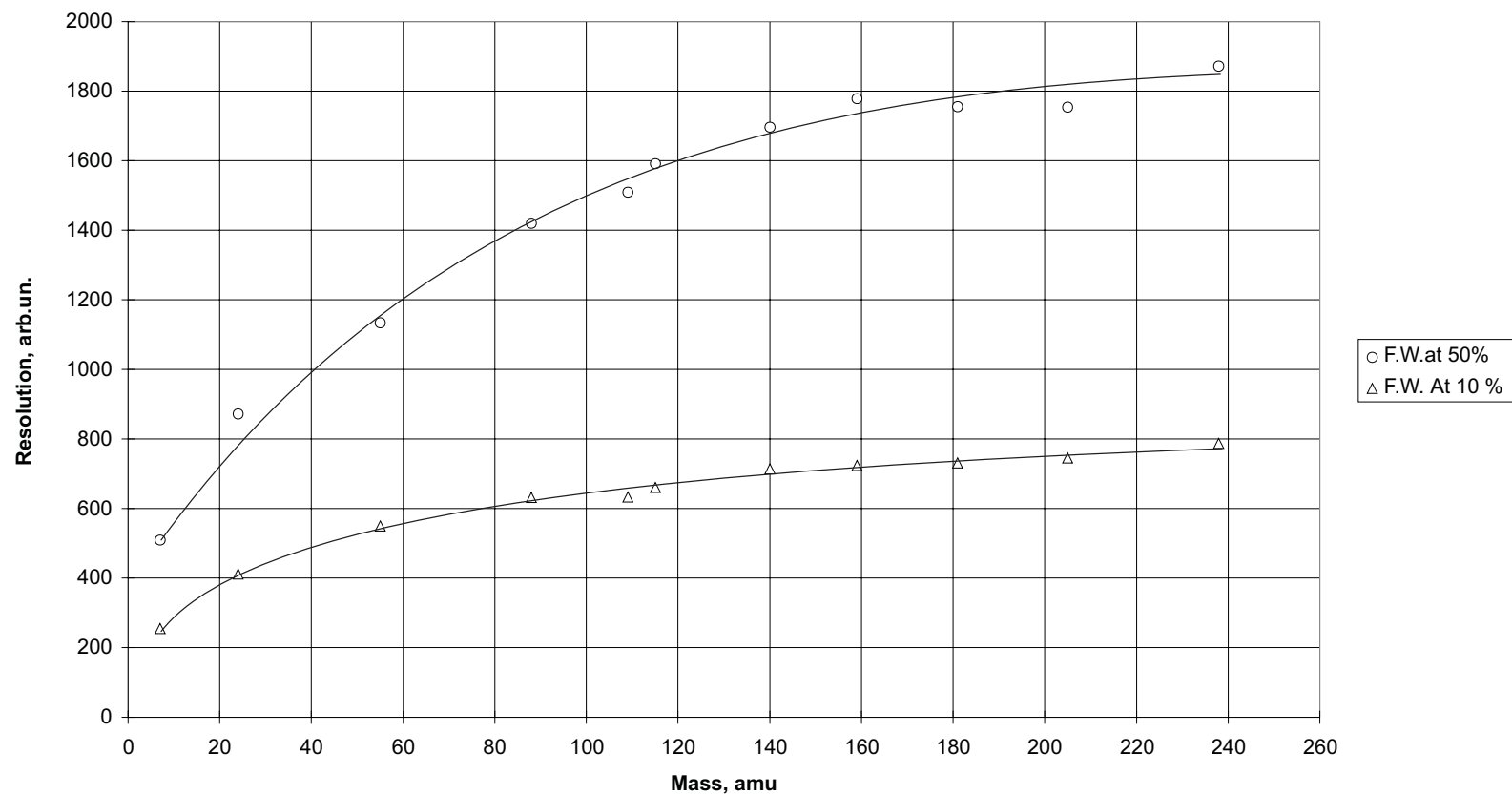

M, amu		7	24	55	88	109	115	140	159	181	205	238
$\Delta M$ , amu	at 10%	0.03	0.06	0.1	0.14	0.17	0.17	0.2	0.22	0.25	0.28	0.3




# 8

## Dynamic range evaluation

- **Dynamic range was evaluated by measuring responses (peak areas) for Ba130, Ba135, Ba137, Ba138 in RO water solutions at 10, 100, 1000 and 15,000 ppb, 3 x 10 s acquisitions (integral of 311,296 OA extractions). At low concentrations Xe130 interference was accounted for (Ba130 = Counts(130)-0.155\*Counts(129))**
- **Correlation factors with linear fit obtained were :**
  - Corr.( TDC) = 0.991**
  - Corr. (TR) = 0.995**

# Linear Dynamic Range

## LINEARITY, Ba

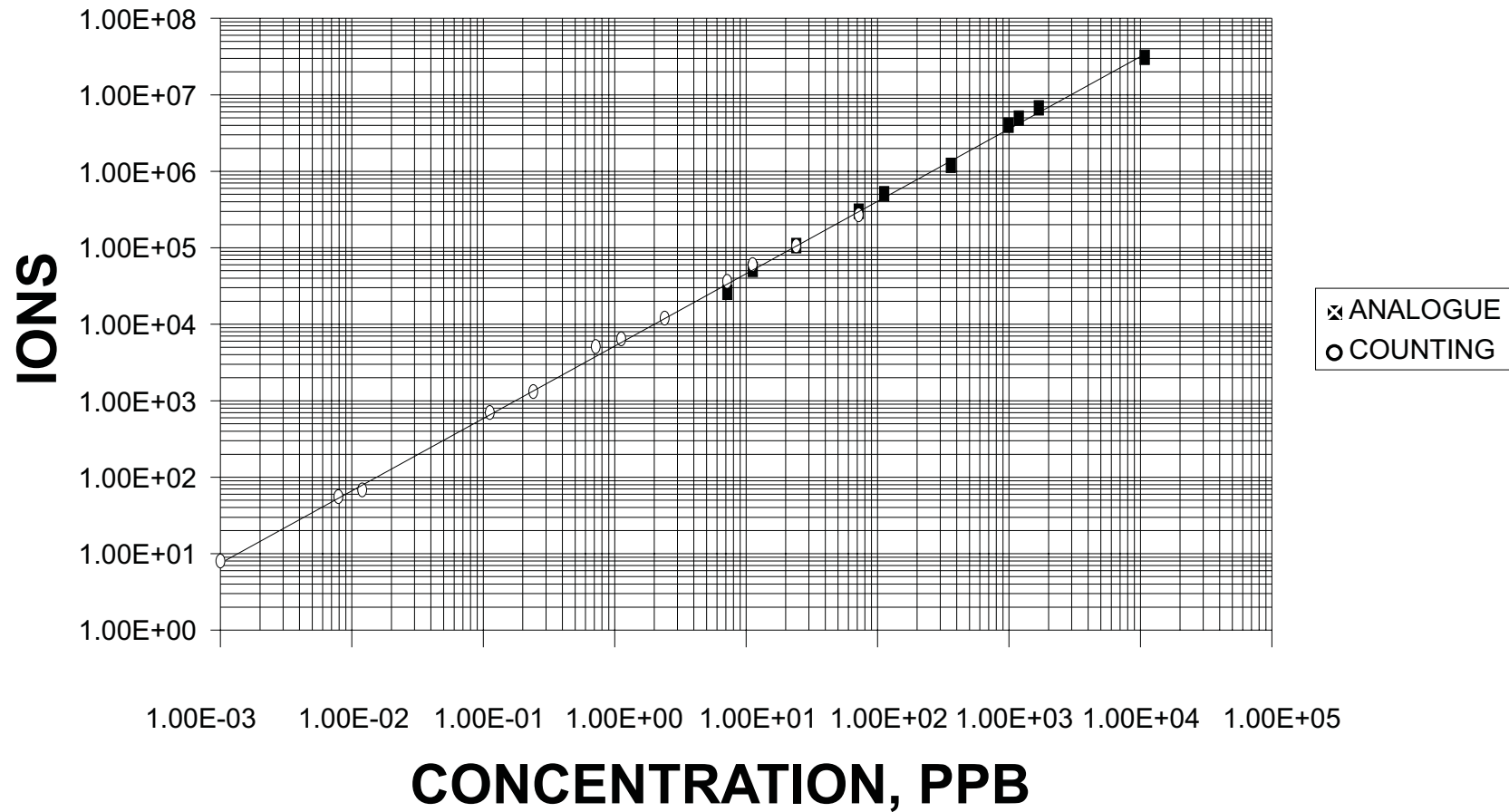

## Detection Limits



**3- $\sigma$  detection limits were determined from the sequence of 10 x 10 s sample-blank acquisitions. Sample: 50 ppt multi-element solution; Blank: RO water blank**

<b>Element</b>	<b>Detection Limit (ng/L)</b>
V, Mn, Co, Rb, Sr, Y, Zr, Nb, Rh, Ag, In, Sc, Ba, Ce, Tb, Ho, Ta, Pb, Bi, U	<1ng/L
Li, Mg, Al, Ti, Cu, Ga, Mo, Pr, Nd, Re, Pt, Au	<10ng/L



**Table 1: Typical DLs for 10 s acquisition with the OptiMass 8000.**

# Detection Limits

Rh at 50 ppt Instrument  $3\sigma$  DL 1 ppt

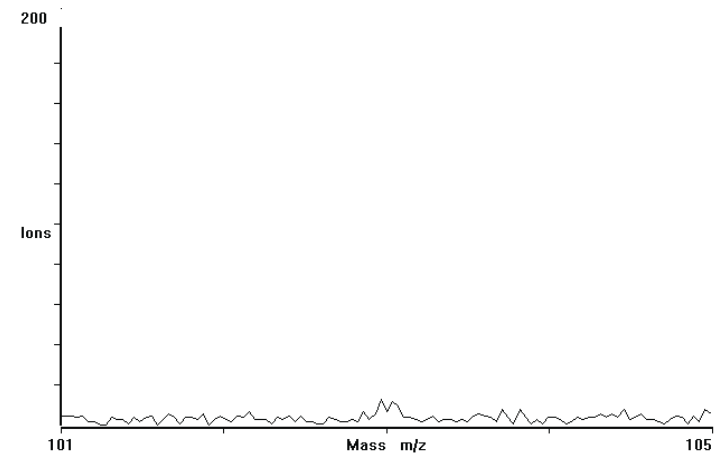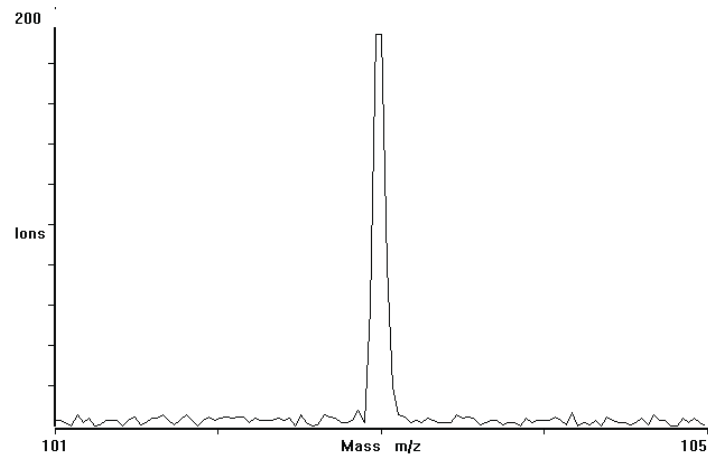

U at 50 ppt Instrument  $3\sigma$  DL 0.5 ppt

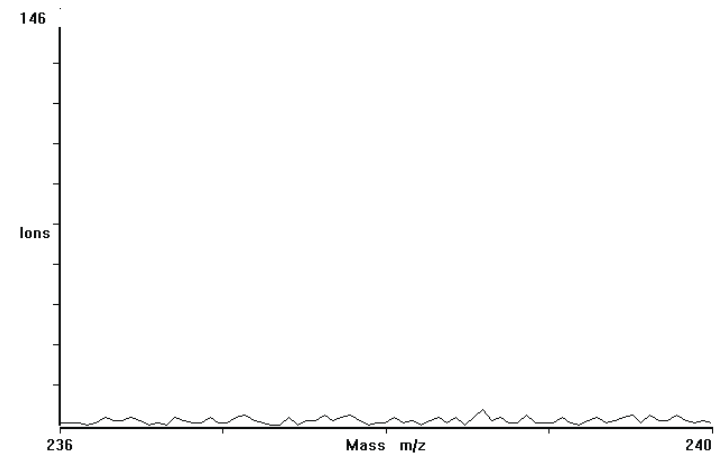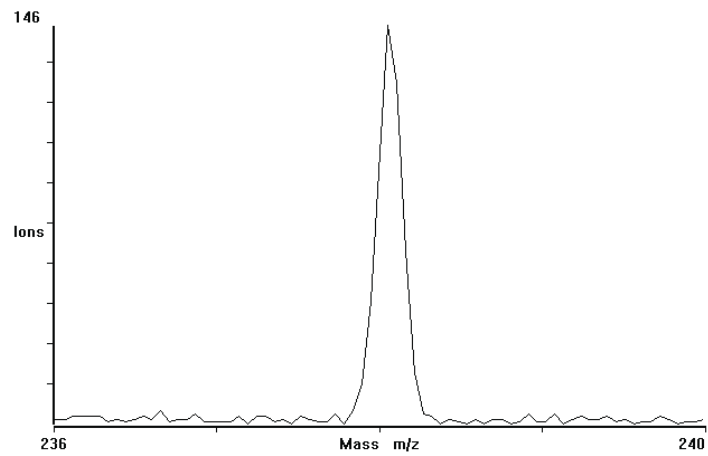


## **Ion Blanking**

**As a time-of-flight analyzer allows all ions into the analyzer, a mechanism is required for preventing highly abundant ions such as  $\text{Ar}^+$  from reaching the detector.**


### **OptiMass 8000 SMARTGATE Ion Blanker**

- effectiveness 100,000 (adjustable)**
- multiple mass blanking rate 10 MHz**
- resolution 25 nS**

## Effectiveness of SMARTGATE Ion Blanker

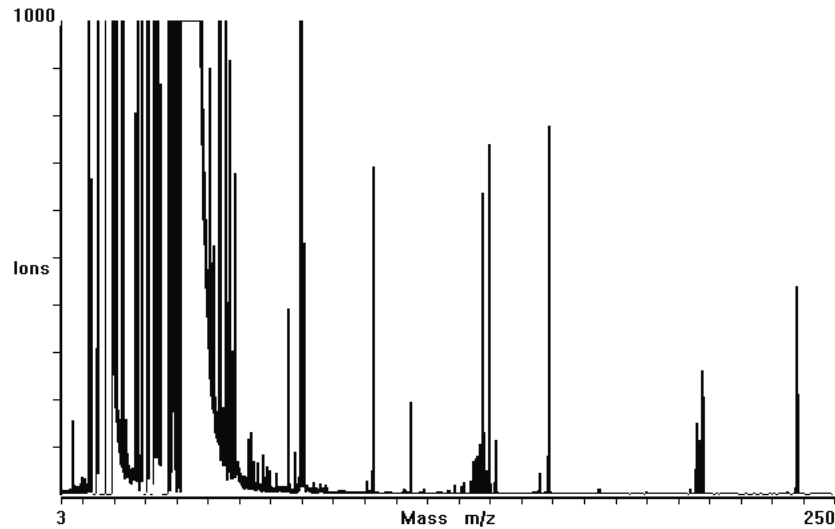


**SMARTGATE: Off**  
**Multiplier voltage: 2.4 kV**

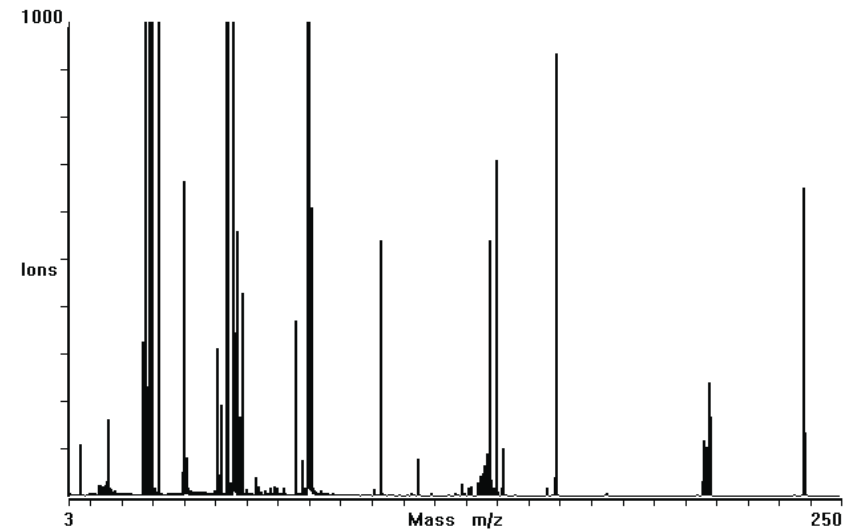

**SMARTGATE: On**  
**Multiplier voltage: 2.4 kV**  
**Ar eliminated to < 2 ppb conc.**

## Evaluating Mass Response Curve

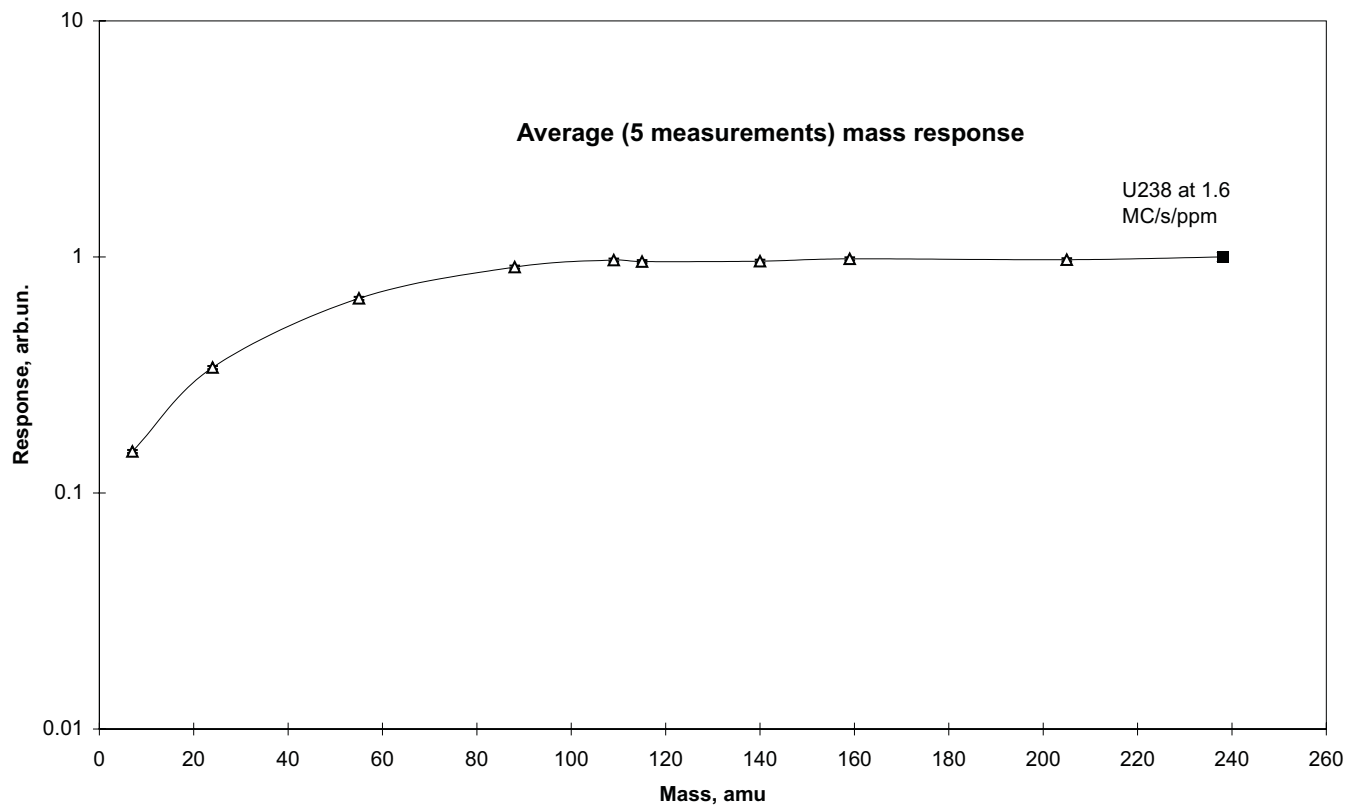

**Mass response curve was measured using 10-element solution at 10 ppb in RO water. Natural abundances and equilibrium degrees of first and second ionization were accounted for.**

15

## **MO/M and M<sup>++</sup>/M<sup>+</sup> Ratios**

**Oxide and double charged ions of Ce and Ba maximal values were defined for 7 x 15 s acquisitions under normal operating conditions.**


**The maximal measured values were :**






16

## CeO/Ce Ratio

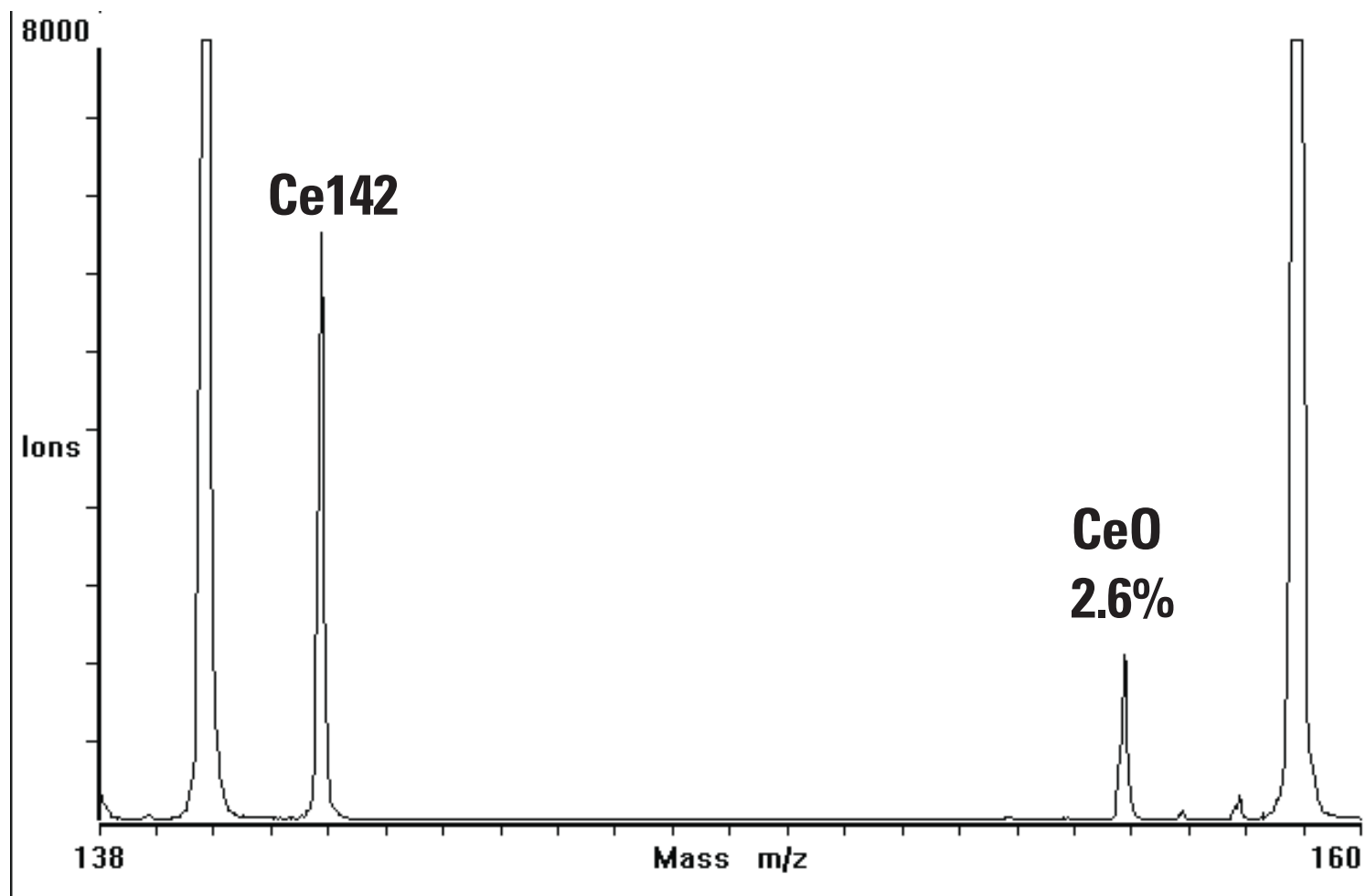

17

# Ba<sup>++</sup>/Ba ratio

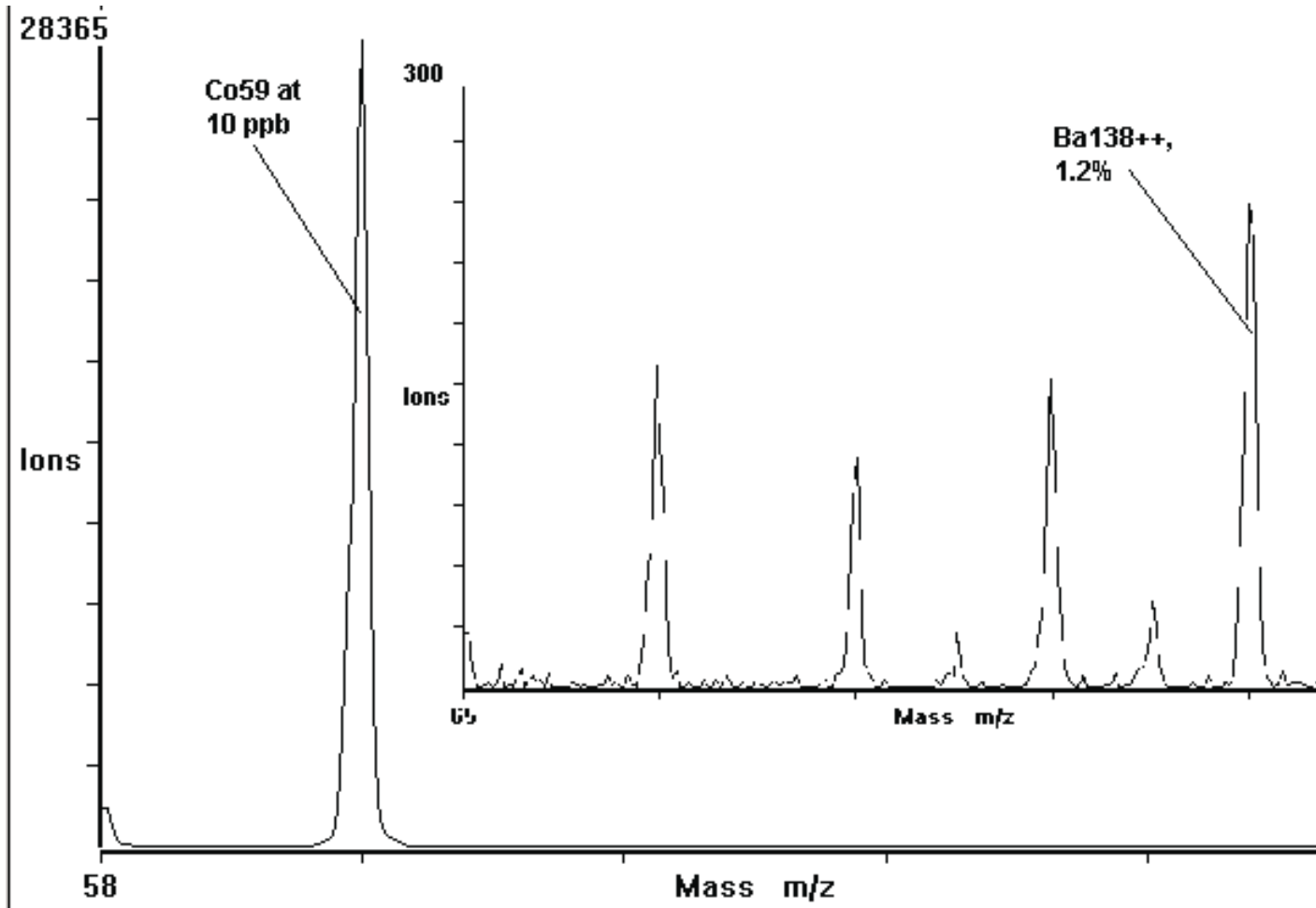

## Ratio Precision

**Ag107/Ag109 ratio precision was evaluated from full mass range measurements using 15 s TDC scans of a 10 ppb solution.**

- **15 s acquisitions:**

- Ag107 — 93,050 counts with 2.3 % RSD;**

- Ag109 — 90,906 counts with 2.2 % RSD;**

- **Integrals for 10 consecutive points:**


- Ag107 — average 931,327 counts at 0.62% RSD;**

- Ag109 — average 911,110 counts at 0.73 % RSD;**

- Ratio RSD — 0.174 %, Statistical limit 0.147% .**

# Ratio Precision

64 x 15 s measurements of 10 ppb Ag solution

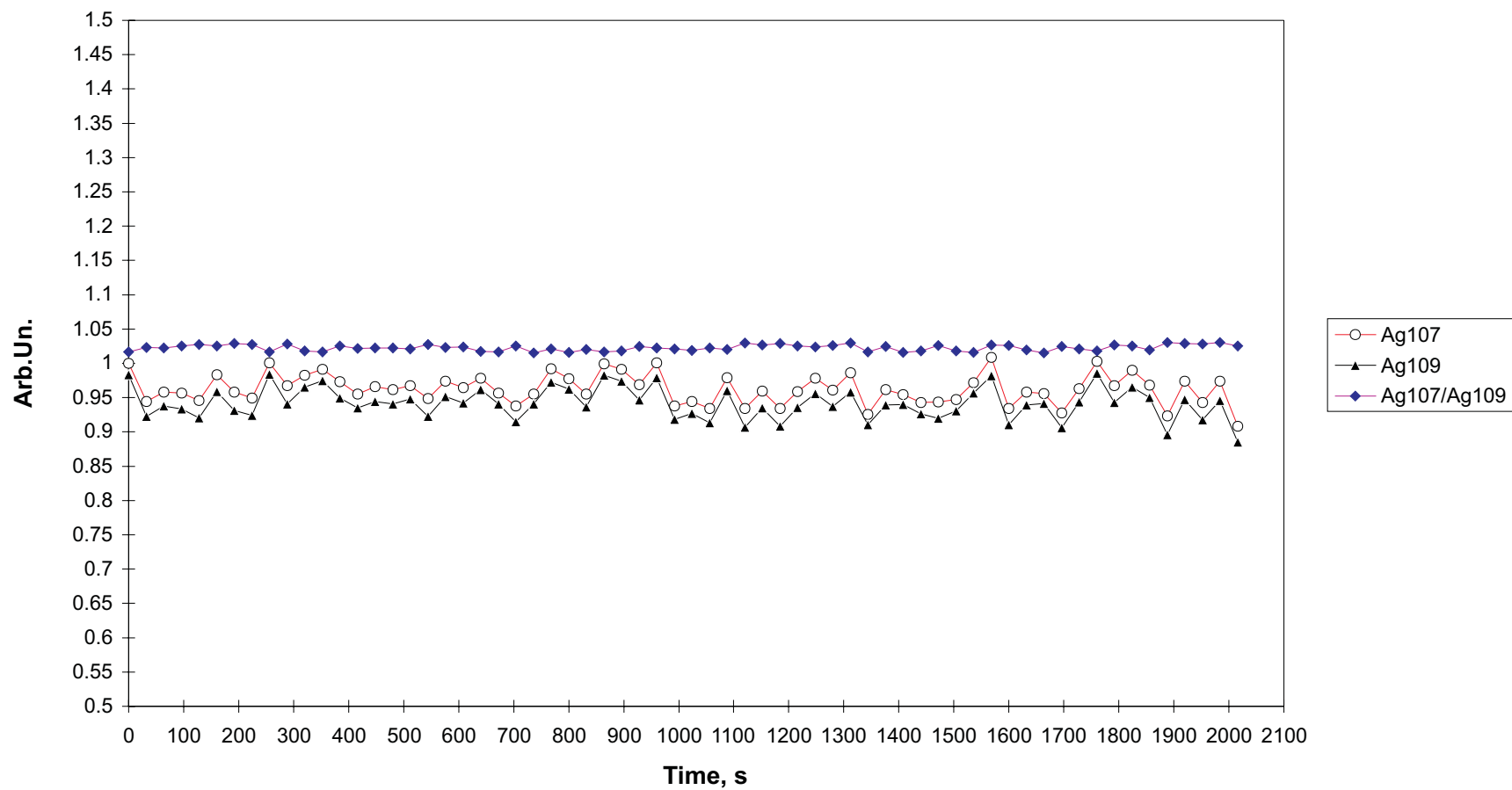

## Conclusions

**OptiMass 8000 under standard operating conditions displays the following performance characteristics:**

- **Resolution  $> 1800$  (f.w.h.m.); 0.03–0.3 a.m.u. At 10 %**
- **Dynamic range  $10^7$**
- **Detection limits 1–10 ppt for 10 s acquisition**
- **Ar elimination efficiency of  $>100,000:1$**
- **Flat mass response for 80–250 a.m.u. Mass range**
- **$M^0/M < 3\%$ ;  $M^{++}/M^+ < 2\%$**
- **Ratio Precision is Approaching Statistical Limit**