

EYELA

SPRAY DRYER MODEL SD-1000

PRINCIPLE

Sample solution is atomized and blown through hot air flow and is collected in the form of dried uniform powder.

ADVANTAGES




Very short drying time enables the treatment of heat sensitive samples.

You can obtain the dried uniform powder within a few minutes.

You can save the time consuming procedure such as concentration or filtration.



TOKYO RIKAKIKAI CO., LTD.
TOKYO JAPAN

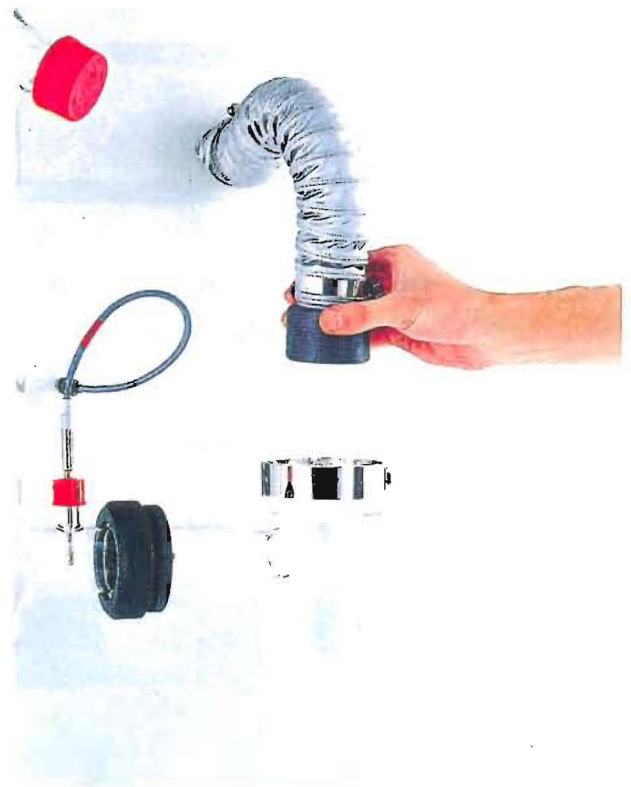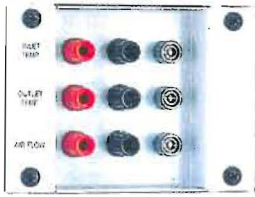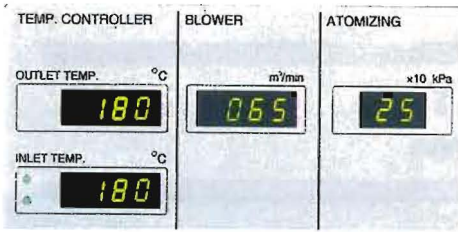


Easy setting/check-up/recording of data


Indicator shows both temp. at inlet and outlet and spraying air pressure. They are also output on recorder

Easy disassemble/cleaning/reassemble

Easily detaches/attaches evaporating tube. One touch coupling of cyclon /suction blower line

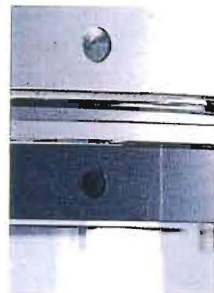

Useful setting mark

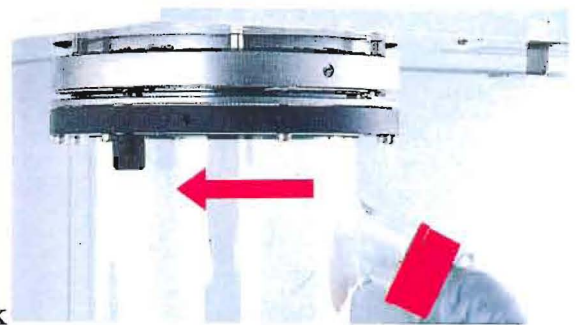

Suction filter at inlet



Airfilter at outlet



Efficient layout of plumbing line


Shortest plumbing line for efficiency.

SPECIFICATIONS

SD-1000

Type	2 way nozzle type Spray nozzle
Drying capacity	Max.approx.1500ml/h
Temp range · accuracy	40~200°C (Temp. at inlet) ±1°C
Dried air volume range	0.2~0.65m ³ /min
Spray air pressure range	49~245kPa (0.5~2.5kg/cm ²)
Feed pump volume range	102~1800ml/h
Sample stirrer rpm · capacity	100~1000rpm · 50ml~2L (water)
Inlet temp. setting · Display	Membrane switch · Digital
Outlet temp.display	Digital
Dried air flow rate setting	Variable resistance · Digital display
Feed pump flow rate setting	Variable resistance
Sprayed air pressure setting	Variable resistance · Digital display
Stirring speed setting	Variable resistance
Safety device	Cover for drying tube, Self-diagnostic device, (Temp. error, Air flow error, Damaged heater, Flow pressure error)
Sprayed air line cleaning device	Automatic by interval timer
Additional function	Sprayed air line valve ON-OFF switch
Recorder output terminal	0~10mV (Inlet temp. · Outlet temp. Dried air flow)
Heater	3kW
Feed pump	Peristaltic pump
Spray nozzle	2 way nozzle (ID 0.4mm)
Air pump for spray	Air compressor for spray (optional)
Drying tube	Hard glass
Cyclone tube	Hard glass
Powder collecting container	Hard glass (600ml)
Stirrer for sample	DC brushless motor 6W, Cobalt magnet
Suction blower	Commutator blower, Max0.65m ³ /min Flow controllable
Dimension of tubing	ID3.15mmxOD5.2mm
Dimension of connector for air for spraying	ID4mmxOD6mm Union for soft urathane
Air pressure for spraying	294kPa (3kg/cm ²), Flow rate more than 25L/min
Dimensions of Exhaust connector	Exhaust OD50mm
Interval timer	OFF or 1~20min (Automatic)
Ambient temp. range for operation	5~35°C
Overall dimensions(mm) · Weight	700Wx620Dx1500H · Approx.110kg
Power requirement	21A, 4.2kVA · AC-200V ø1, 50/60Hz

FLOW DIAGRAM

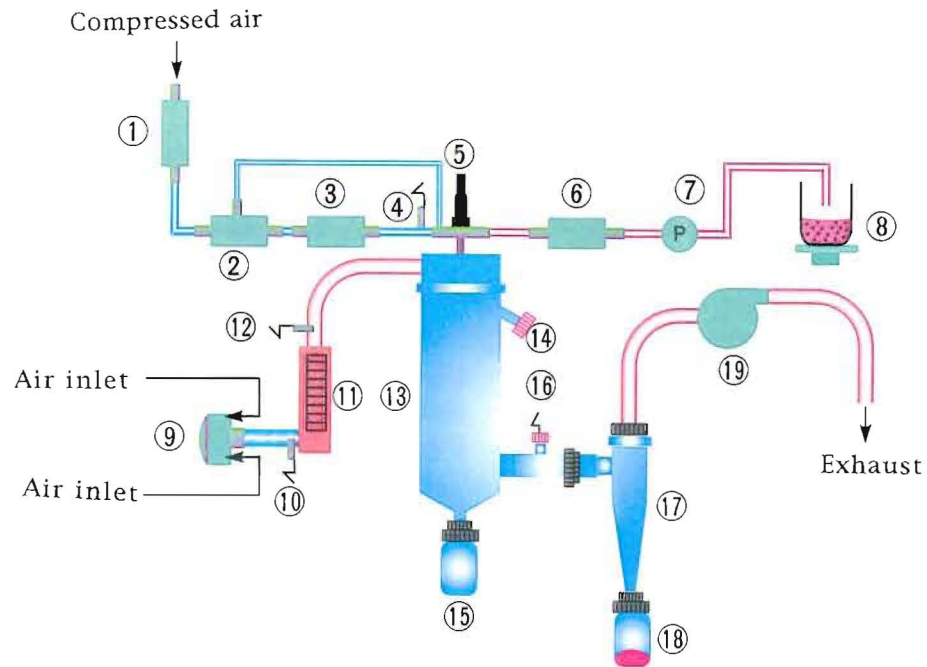

1. Sprayed air line valve (Solenoid valve)
2. 3 way solenoid valve
3. Regulator
4. Pressure sensor (sprayed air)
5. Spray nozzle
6. Pump pressure sensor
7. Pump
8. Sample stirrer (magnetic stirrer)
9. Suction filter
10. Sensor (wind)
11. Heater
12. Sensor (Inlet temp.)
13. Drying tube
14. Spray nozzle purification orifice
15. Separator
16. Sensor (outlet)
17. Cyclone tube
18. Powder collecting flask
19. Suction blower

TOKYO RIKAKIKAI CO.,LTD.

Nihonbashi Hon-cho Bldg. 6F, 3-3-4 Hon-cho, Nihonbashi, Chuo-ku, Tokyo, Japan

Ph: 81/3-5201-6462 Fax: 81/3-3241-0177